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Zafer Bayram

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Shimadzu Corporation's new ultra high-performance liquid chromatography (UHPLC) system, Nexera, delivers high-speed and high-resolution analysis without sacrificing basic performance parameters such as excellent precision (+/-006 % RSD), near-zero carryover (0,0015% max), good linearity, stability and repeatability (0,25 % RSD max). Offering these attributes enables Nexera to provide superior performance in various fields, covering high-resolution and high-speed separation, as well as high-sensitivity detection.

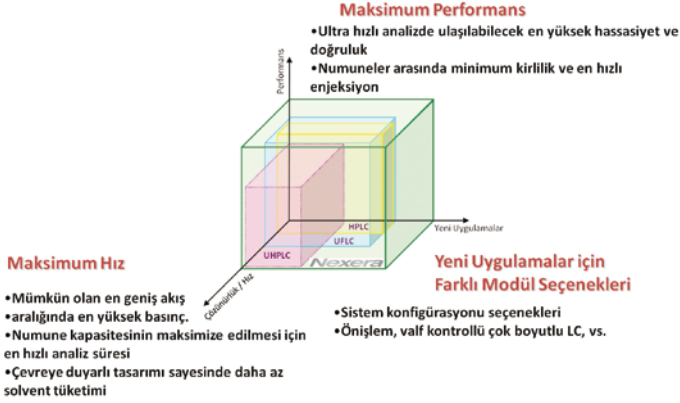
Development Concept Of Nexera: To Produce Next Generation UHPLC Without Compromise Of Performance

Achieving new, greater performance levels, the next-generation Nexera LC from Shimadzu addresses all analysis requirements: conventional, high speed-ultra-high speed, high resolution or high temperature, with no compromise in data integrity.

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
Development Concept of Nexera

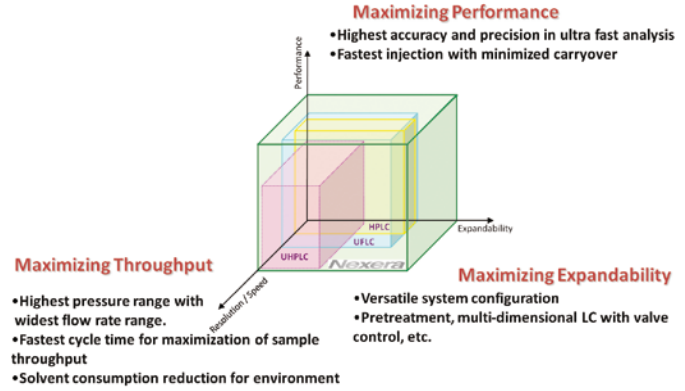To produce next generation UHPLC without compromise of performance

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
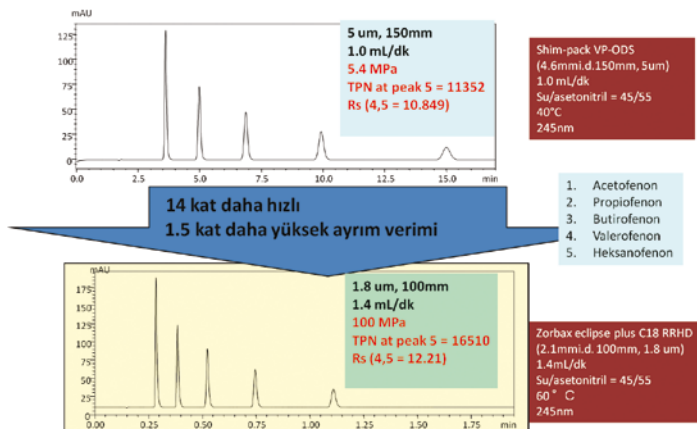
Speed and Resolution

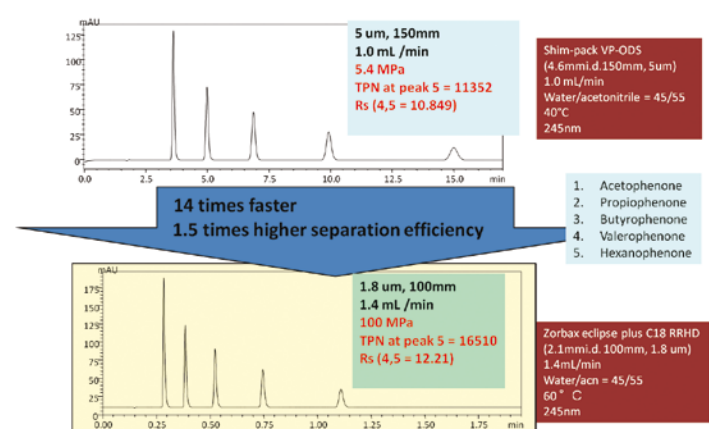Maximizing Performance and Throughput

Nexera offers precise solvent delivery, excellent reproducibility and near-zero carryover, and ensures fully reliable analysis in all application fields through refinement of each component.

Thanks to its S4 Plunger (Super Smooth Surface Sapphire), high strength seal material and new high pressure valve design; Shimadzu's Nexera is optimized to enable analysis at pressures up to 130 MPa (18.850 psi). Without compromising performance and data quality, the maximized pressure over a wide flow range enables various types of analyses, including micro LC to conventional LC, regular pressure LC to ultra high-pressure LC, and isocratic LC to ternary high-pressure gradient and multi-dimensional LC. The system is also suitable for LCMS and LCMSMS applications.

Nexera achieves a perfect balance between the highest resolution and highest speed. Highest resolution in a very short interval is accomplished by Nexera's ultra-high pressure resistance for a broad flow range and the world's fastest sample injector. Ultra-high pressure (up to 130 MPa) fluidic modules can utilize all the high-resolution column. Nexera facilitates the use of long small-particle columns to analyze trace-level compounds in complex samples.

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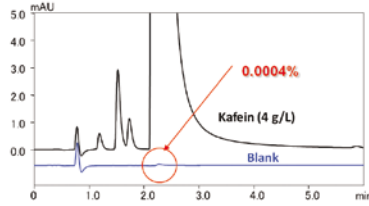
Carryover - Caffeine -


Carryover of neutral compounds (Caffeine) @ 100 MPa

Specification : < 0.0015%

Results : 0.0004 %

Without injection port rinse, needle internal rinse and needle outer surface rinse


	Carryover
Caffeine 2000 ng/µL	—
Post-blank 1	0.0004%
Post-blank 2	N.D.

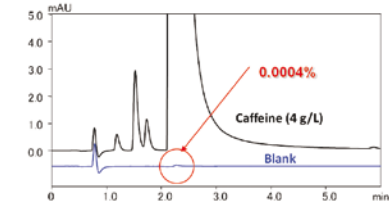Column : ODS (2.0 mmI.D. x 100 mmL, 1.8 µm)
 Mobile Phase : Methanol, Water= 2/8
 Flow rate : 0.4 mL/min
 Temperature : 40 °C
 Injection volume : 5 µL

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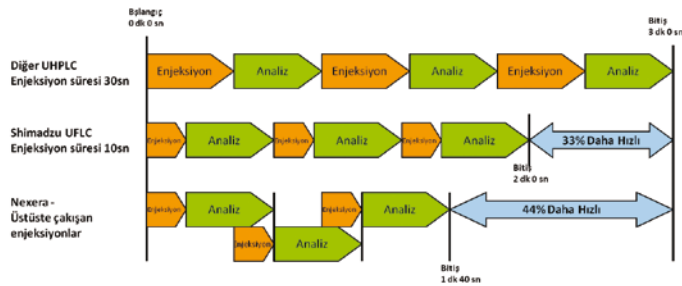
Shimadzu continues to improve its low-carryover reputation with the Nexera autosampler. With reduction of the needle contact area, special coatings, surface treatments and a new needle seal, the SIL-30AC autosampler reaches a new level of low carryover performance.

In addition to minimizing the fundamental causes for compound adsorption, Nexera offers multiple rinse modes with multiple rinse solvents to eliminate carryover of even the most stubborn of compounds. Active rinsing of all contact surfaces can be selected.

The autosampler also offers fixed-loop injection for ultra-fast separation with minimized peak width and gradient delay. An overlapping injection function is standard with Nexera to allow for a maximum reduction in cycle time.

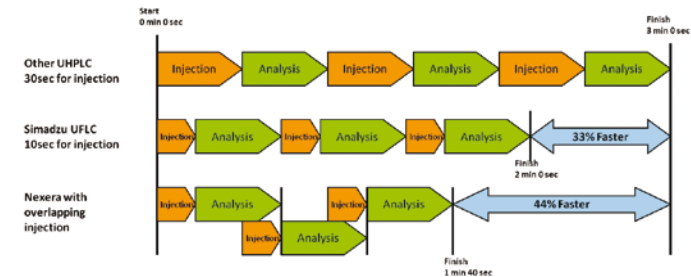Overlap Injections

For high-temperature LC analysis, Nexera offers efficient pre-heating for solvents to enable stable and reliable conditions. The advanced Intelligent Heat Balancer (IHB) minimizes band broadening during high-temperature analysis through the use of independently controlled heat zones within the column compartment, ensuring accurate column temperature regardless of flow rate. The post-column cooler can reduce detector noise when equipped for high-temperature analysis up to 150°C with a temperature setting range of +/-0,05°C.

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
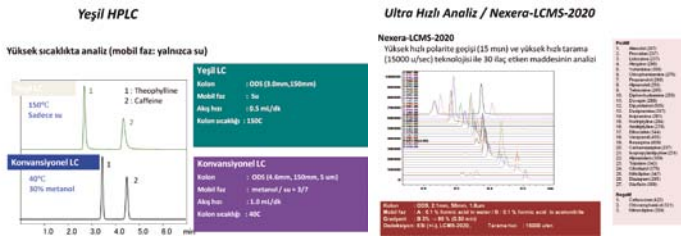
Maximizing Expandability

The Nexera column oven and autosampler, in concert with the modular nature of the system, expand the application range. These applications include high-temperature analysis, green LC, auto-sample pretreatment and multidimensional LC separation.

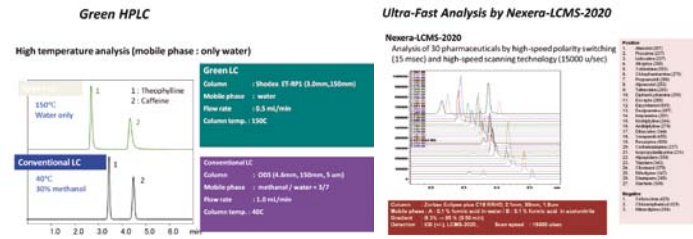Each Nexera component has undergone core improvements to achieve unmatched system performance. In an era when a higher standard of performance is expected, Nexera responds to our customers' needs in all analysis fields: conventional, high speed - ultra-high speed, high resolution, high temperature, green LC, multi-dimensional LC. The UHPLC system that meets the needs of a wide range of applications is none other than Nexera.

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"At Shimadzu, our passion has always been to improve performance, ease of use and flexibility for all HPLC environments," said Shuzo Maruyama, General Manager of LC Business Unit. "That passion continues with Nexera, the only no compromise UHPLC system to offer the reliability, versatility, and superior basic performance for all applications."

About Ant Teknik

Ant Teknik, founded in 1999 in Istanbul, is offering analytical instruments, technical service and spare parts; validation, application and training as well turn-key laboratory project services for Quality Control and Research&Development laboratories. Today, we offer innovative analytical solutions for various fields such as pharmaceutical, food, chemistry and environment. We have offices in Istanbul, Ankara and Izmir; about 60 employees; and resellers servicing to 15 different regions.

We represent the most respected producers of laboratory instruments like Shimadzu Corporation, AB Sciex, Celsis and JeioTech in Turkey; while we provide chromatography consumables from well-known suppliers like GLSciences, Kromasil and PSS.

As Ant Teknik family, we continue to share our experiences with you to contribute to your success.